

## **SAFETY FEATURES**

**To prevent entrapment the buddy-system must be used.**

**Two individuals must be present each time the bulk sterilizer is started.**

**The inside of the autoclave is observed to confirm that no one is inside prior to activating.**

**Both individuals will remain outside the autoclave to ensure that no one is inside as the door is closed and the sterilizer cycle is being activated.**

**By design the door to the sterilizer only closes very slowly to reduce the possibility of entrapment.**

**Emergency Shut-Off (red button) located on both sides of the autoclave.**

**Emergency Electrical Cut-Off located to left of unit on prep-side.**

SC #259  
Eff 7/19